

# Microscope Decontamination Guide

## Personal Protective Equipment

- Gloves, goggles, lab coat are optional for BL1
- Please use **FRESH GLOVES**, located by sinks
- Notify us of vinyl allergy
- BL2 requires additional PPE and Training**


## Upon Completion

- Clean oil from lenses
- Wipe-down as instructed
- Wipe and dry eyepieces, keyboard, mouse
- Turn off scope using 70% ethanol wipe

## Equipment Wipe-Down


### ① Decontaminate

- Spray large kimwipe with 70% ethanol
- Wipe stage and touch-points on microscope including joystick
- Wipe eyepiece and its pillar, computer keyboard, and mouse
- Use ethanol kimwipe to **TURN OFF** power switches and laser keys
- Wipe **EVEN IF** you used gloves

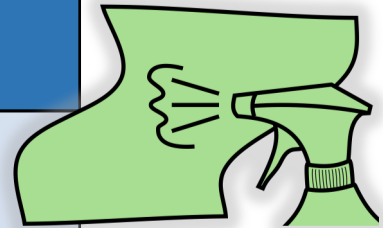

### ② Dry

- Use fresh kimwipe
- Dry eyepieces, scope body, keyboard, and mouse.

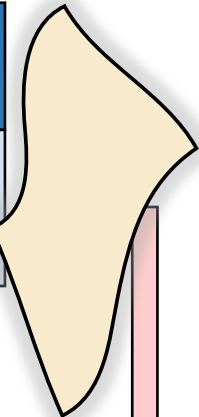

### ③ Dispose

- Collect kimwipes
- Dispose of gloves and kimwipes in biosafety box

